## **Instruction Manual Manuel d'instructions Manual de instrucciones Bedienungsanleitung**

## **Fisherbrand Infrared Vortex Stirrer 3000 t/min Part Number: 1328-4769**

### **General Information / Informations Générales / Información General / Allgemeine Hinweise**

Before using the unit, please read the following instruction manual carefully.

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.

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#### **This unit must be used for laboratory applications only.**

The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions.

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#### **This unit has been designed and manufactured in compliance with the following standards:**

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Safety requirements for electrical equipment for measurement, control and for laboratory use

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**IEC/EN 61010-1  
IEC/EN61010-2-051**

Electrical equipment for laboratory use

General requirement - Canadian electrical code

**UL 61010-1  
CAN/CSA-C22.2 No.61010-1**

Fisher Scientific reserves the right to modify the characteristics of its products with the aim to constantly improving their quality.

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The plug disconnects the instrument. Therefore, place the instrument where it can be quickly disconnected.

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The heated solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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Position the instrument on a flat surface, with a distance from the wall of 30 cm (at least).

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it.

It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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# Contents / Index / Índice / Inhalt

## EN

1.	INTRODUCTION.....	4
2.	ASSEMBLY AND INSTALLATION.....	4
	2.1 ELECTRICAL CONNECTIONS.....	4
	2.2 START-UP.....	4
	2.3 INFORMATION ABOUT CONSTRUCTION MATERIALS.....	4
3.	OPERATING CONTROLS.....	4
4.	MAINTENANCE.....	5
	4.1 CLEANING.....	5
5.	TECHNICAL DATA.....	5
6.	ACCESSORIES.....	5

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Mixing by spinning can be carried out quickly and safely by placing the tube containing the sample on the mixing cup head. The Vortex Mixer can be used in automatic (TOUCH) or continuous mode (CONT.): in both cases the mixing speed can be adjusted using the speed selector knob (1).

When the instrument is switched on, it is always set in TOUCH mode. Mixing starts when the test-tube crosses the infrared sensor (6) field near the mixing cup head (4). For continuous mode press the mode select button (2), mixing starts automatically at the selected speed, no pressure is required on the cup head.

The led (5) indicates that the instrument is running in TOUCH mode.

The press-on fitting means that the mixing cup head can easily be replaced by any of the various accessories available for test-tubes of different sizes.

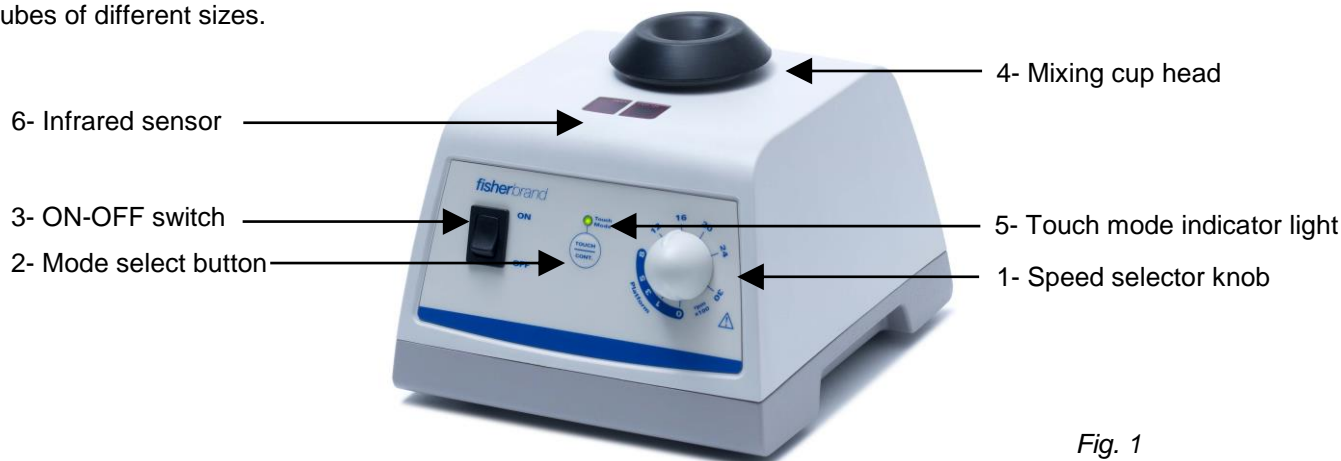
Fig. 1

## 2. Assembly and installation

Check the integrity of the unit after unpacking. The box includes:

- Fisherbrand Infrared Vortex Stirrer 3000 t/min with mixing cup head
- EU and UK adapter plug
- Power supply 100-240V/12V no plug
- Instruction manual

### 2.1 Electrical connections

After having unpacked the instrument, place the unit on the laboratory bench.

Before connecting the instrument to the power supply, make sure that the values on the rating plate correspond to those of the power supply. Connect the unit to the power supply using the transformer supplied.

Ensure that the socket and the relative cut-off device conform to current safety norms and are easy to reach.

### 2.2 Start-up

Check that the mixing cup head is correctly positioned. Also check that the speed regulation knob is turned to "0 rpm". Use the on-off switch on the front of the unit to turn the unit on (position "ON"). The start-up and the control of the speed are effected with the knob. When switched on the unit starts up in "TOUCH" mode with the indicator lights on.

### 2.3 Information about construction materials

Housing	Zinc alloy/Polymer	Feet	PVC
Paint	PET	Platform	PP
Frontal control panel	PET	Foamed part for platforms	PE
Cup head	SEBS		

## 3. Operating controls

Connect the unit to mains and turn it on using the on-off switch. Select the operating mode using the mode select button on the front of the unit.

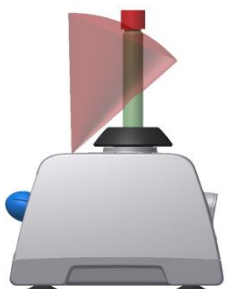
Fig. 2

When the **Touch** mode is selected, mixing starts automatically when the test-tube crosses the sensor field (infrared optical system) shown in Figure 2.

In **Continuous** mode the mixing cup head runs continuously at the set speed.

In both operating modes there is no need for any pressure on the mixing cup head. The sensor field is purpose-designed to detect the presence of any kind of test-tube. Vortex Mixers carry-out mixing by spinning without exerting pressure on the cup head, thus reducing effort by the operator. However, it is important to grip the tube firmly during mixing.

## CONTINUOUS-TOUCH

Pushing the mode select button, the respective modes are activated:

- **Touch** (indicator light on) allows automatic operation mode even with different accessories.
- **Continuous** (indicator light off). The instrument runs in continuous mode with various accessories.

## SPEED REGULATION KNOB

The knob on the front of the unit can be used for quick precision regulation of mixing speed from 0 to 3000 rpm.

## ON-OFF SWITCH

The on-off switch turns the unit on and off. If the switch is in the "OFF" position the unit is off; if the switch is in the "ON" position the unit is on.

Always turn the unit off after use.

## 4. Maintenance

No routine or extraordinary maintenance is necessary apart from periodically cleaning the unit as described in this manual. In compliance with the product guarantee law, repairs to our units must be carried out in our factory, unless previously agreed otherwise with local distributors.

To replace the cup head, pull it upwards and off. The other push-fit accessories can then be inserted.

### 4.1 Cleaning

Disconnect the unit from the power supply and use a cloth dampened with an non-inflammable non-aggressive detergent.

## 5. Technical data

Power supply input	AC 100-240V ( $\pm 10\%$ ); 50-60 Hz ; 0.5A
Power supply output	DC 12V ; 1.25A
Dimensions (WxHxD)	150x130x165 mm (5.9x5.1x6.5 in)
Weight	2.7 Kg (6.2 lb)
Speed range	0-3000 rpm
Spin diameter	4.5 mm (0.177 in)
Max. weight on the mixing cup head / Accessories	0.5 kg (1.1 lb)
Permitted operation	Continuous
Operation mode	Continuous or Sensor
Level of electrical protection IEC/EN60529+A1	IP 42
Noise level	<< 80 dBA
Admitted temperature	Storage: -10...+60 °C (+14...+140 F°) Working: 5°...+40 °C (+41...+104 °F)
Admitted humidity	Max. 80%
Power	15 W
Overvoltage category	II
Pollution degree IEC/EN61010-1	2
Max. altitude	2000 m

## 6. Accessories

13256959	Foam stand for 19 microvials 1.5 ml - Eppendorf®
11792573	Customizable soft foam top
11712583	Foam stand 5 test tubes Ø16mm
11792583	Foam stand for microtiter
11702593	Small rubber supporting plate Ø 50mm
11840992	Foam stand 4 test tubes Ø29mm

		
13256959	11792573	11712583
		
11792583	11702593	11840992

Do not exceed a speed of 800 rpm when any one of the above accessories is in use. High speeds may cause the sample to spill.

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